

### RFID Sticker for Improved Traceability



Smart Tags are typically affixed to pipettes or beakers to track the sample more easily.

#### **Choosing an area to fix the Smart Tag**

The area must be:

- flat (if only round areas are available, choose a place with a large radius)
- far from metal
- away from the gripping area
- easy to hold next to the RFID reader

#### **Fixing the Smart Tag to the container**

- 1 Clean the area with a mild solvent, e.g., isopropanol or methanol.
- 2 Wait until the area is completely dry.
- 3 Fix the Smart Tag to the area.

**Mettler-Toledo GmbH**

Im Langacher 44  
8606 Greifensee, Switzerland  
[www.mt.com/contact](http://www.mt.com/contact)

Subject to technical changes.  
© Mettler-Toledo GmbH 10/2019  
30104632C en



30104632